

Roti®-DipSlides

Sterile dip slides with solid, ready-made culture medium for easy microbiological analysis of solutions and surfaces.



ready-to-use

S



Tried-and-tested media in a new look –
Now with optimized plastic carriers!

Advantages

- 2 in 1: DipSlide and contact agar
- Very easy application
- Reliable assay and easy interpretation
- Detection limit is 10² cfu/ml solution
- Acc. to ISO 11133
- Detailed instructions-for-use included

Characteristics


- Made of a flexible plastic carriers, sterile
- Coated with solid culture medium on both sides
- Total analysis area is 11,44 cm² per side
- After dipping in the solution or pressing on the surface, the slide is placed back into the sterile tube and incubated
- Storage temp.: +4 to 15 °C

Suitable for:


- Microbiological laboratories
- Hygiene control in food production and sports sector
- Routine checks
- Inspection of air conditioning

Application


- ▶ Remove carrier from sterile vial. The screw cap automatically serves as a holder
- ▶ Press agar gently onto surface or immerse in liquid
- ▶ Insert the carrier back into the tube and screw it
- ▶ Incubate at a temperature suitable
- ▶ Evaluate

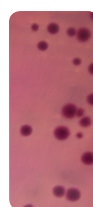


VRB

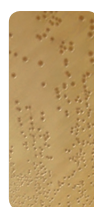

PC

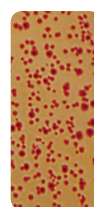

TTC-Agar

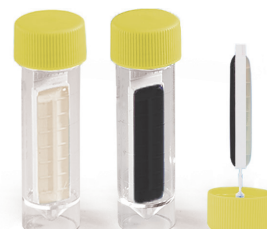

For easy interpretation of results, use the following table

No. of colonies	Result	Interpretation
0	Negative	No contamination
1–10	Weakly positive	Weak contamination
11–20	Positive	Low contamination
21–100	Strongly positive	High contamination
>100	Highly positive	Very high contamination

Important information on refrigerated transports

All products identified as being particularly temperature-sensitive are shipped in **special ice Boxes with freezer packs or in dry ice.**

Additional costs resulting from such, **will be invoiced** (further information on request).

Please note:

In order to guarantee optimum product quality, the dispatch of refrigerated transport products will be carried out **on Mon and Tue only outside Germany!** Slight delays in delivery are therefore possible.

Roti®-DipSlides

ready-to-use

S



ROTI®-DipSlide MacCon/Cetr

For detection of enterobacteria like salmonellae or shigellae and pseudomonads in liquids or on surfaces.

Art. No.	Pack Qty.	Packaging
8045.1	20 pieces	box

ROTI®-DipSlide Nutrient-TTC

For total germ count in liquids or on surfaces.

Art. No.	Pack Qty.	Packaging
2837.1	20 pieces	box

ROTI®-DipSlide CASO-TTC/RBCenr

For total germ count and for detection of yeasts and moulds in solution or on surface.

Art. No.	Pack Qty.	Packaging
3934.1	20 pieces	box

Suitable for analysis of water-mixable cooling lubricants, air conditions, water-soluble paint, and cosmetics.

ROTI®-DipSlide PCA/VRB

For total germ count and for detection of Coliforms in liquids or on surfaces.

Art. No.	Pack Qty.	Packaging
2925.1	20 pieces	box

ROTI®-DipSlide PCA

For total germ count in liquids or on surfaces.

Art. No.	Pack Qty.	Packaging
2927.1	20 pieces	box

ROTI®-DipSlide PCA/Legionella

For total germ count and for detection of legionells in solution or on surface.

Art. No.	Pack Qty.	Packaging
2493.1	20 pieces	box

Suitable for supervision of units for heating of drinking water (Drinking Water Regulations), for surveillance of ventilation and air conditioning, as well as for monitoring of the microbiological status of evaporative cooling systems.

ROTI®-DipSlide PCA/RBCplus

For total germ count and for detection of yeasts and moulds in liquids or on surfaces.

Art. No.	Pack Qty.	Packaging
2926.1	20 pieces	box

ROTI®-DipSlide D/Enneutral

For total germ count on surface after disinfection. For hygiene monitoring. Contains 5 neutralizers.

Art. No.	Pack Qty.	Packaging
8023.1	20 pieces	box

ROTI®-DipSlide Sab4plus

For detection of yeasts and moulds in liquids or on surfaces.

Art. No.	Pack Qty.	Packaging
2956.1	20 pieces	box

ROTI®-DipSlide PCA/TSA-Letheen

For total germ count in liquids or on surfaces. Suitable for analysis of cosmetics and of preservative containing products or after disinfection. Contains two neutralizers.

Warning H319

Art. No.	Pack Qty.	Packaging
8715.1	20 pieces	box

ROTI®-DipSlide PCA/VRBD

For total germ count and for detection of enterobacteria in liquids or on surfaces.

Art. No.	Pack Qty.	Packaging
2924.1	20 pieces	box

ROTI®-DipSlide CLED/MacCon

For detection of bacteria of the urinary tract and of enterobacteria/Coliforms in liquids or on surfaces.

Art. No.	Pack Qty.	Packaging
2957.1	20 pieces	box

ROTI®-DipSlide VRBD

For detection of enterobacteria in liquids or on surfaces.

Art. No.	Pack Qty.	Packaging
2954.1	20 pieces	box

Kontakt Deutschland:

Bestellungen zum NULLTARIF 0800 5699-000
Tel.: 0721 5606-0 · Fax: 0721 5606-149
bestellungen@carlroth.de · www.carlroth.de

Carl Roth GmbH + Co. KG
Schoemperlenstr. 3-5 · 76185 Karlsruhe

Es gelten die allgemeinen Geschäftsbedingungen der Carl Roth GmbH + Co. KG.

Contact international:

Phone +49 721/5606 510
Fax +49 721/5606 111
www.carlroth.com

Carl Roth GmbH + Co. KG
Schoemperlenstr. 3-5 · 76185 Karlsruhe

Export prices might be higher. All supplies and deliveries are subject to the terms and conditions of sale and delivery of Carl Roth GmbH + Co. KG, Karlsruhe.