


# ROTI® DipSlides

Sterile dip slides with solid, ready-made culture medium for easy microbiological analysis of solutions and surfaces.

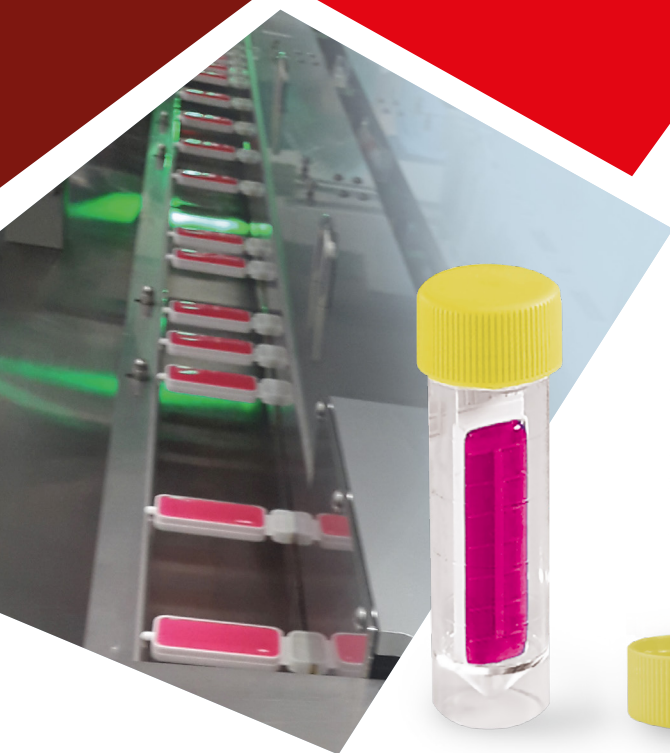

**S** ready-to-use

Tried-and-tested media in a new look –  
Now with optimized plastic carriers!

#### Advantages

- **2 in 1:** Immersion culture media and impression agar
- Easy to use
- Reliable assay and easy interpretation
- Suitable for determining the microbiological contamination and for hygiene controls in a variety of domains
- Including detailed instructions-for-use

#### Characteristics


- Made of a flexible plastic carriers, sterile
- Coated with solid culture medium on both sides
- Total analysis area is 11,44 cm<sup>2</sup> per side
- After dipping in the solution or pressing on the surface, the slide is placed back into the sterile tube and incubated
- Storage temp.: +15 bis +25 °C

#### Application

- Remove carrier from sterile vial. The screw cap automatically serves as a holder.
- Press agar gently onto surface or immerse in liquid
- Insert the carrier back into the tube and screw it
- Incubate at a temperature suitable
- Evaluate

#### Suitable for:

- Microbiological laboratories
- Hygiene control in food production and sports sector
- Routine checks
- Inspection of air conditioning

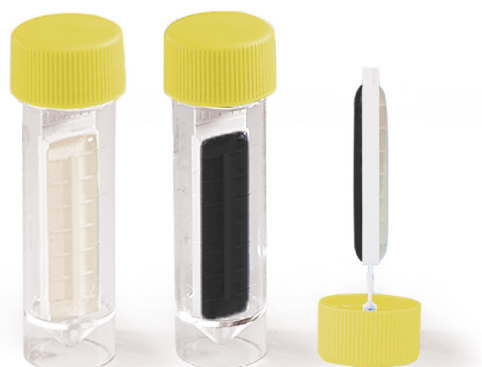

**For easy interpretation of results,  
use the following table:**

No. of colonies	Result	Interpretation
0	Negative	No contamination
1–10	Weakly positive	Weak contamination
11–20	Positive	Low contamination
21–100	Strongly positive	High contamination
>100	Highly positive	Very high contamination

**Detection limit is 10<sup>2</sup> cfu/ml solution assayed.**  
For detailed instructions on interpretation see instruction-for-use.

# ROTI®DipSlides

**S****ready-to-use**

## ROTI®DipSlide MacCon/Cetr

For detection of enterobacteria like salmonellae or shigellae and pseudomonads in liquids or on surfaces.

Art. No.	Pack Qty.	Pack.
8045.1	20 unit(s)	box

## ROTI®DipSlide Nutrient-TTC

For total germ count in liquids or on surfaces.

Art. No.	Pack Qty.	Pack.
2837.1	20 unit(s)	box

## ROTI®DipSlide CASO-TTC/RBCenr

For total germ count and for detection of yeasts and moulds in solution or on surface. Suitable for analysis of water-mixable cooling lubricants, air conditions, water-soluble paint, and cosmetics.

Art. No.	Pack Qty.	Pack.
3934.1	20 unit(s)	box

## ROTI®DipSlide PCA/RBCplus

For total germ count and for detection of yeasts and moulds in liquids or on surfaces.

Art. No.	Pack Qty.	Pack.
2926.1	20 unit(s)	box

## ROTI®DipSlide D/Eneutral

For total germ count on surface after disinfection. For hygiene monitoring. Contains 5 neutralizers.

Art. No.	Pack Qty.	Pack.
8023.1	20 unit(s)	box

## ROTI®DipSlide Sab4plus

For detection of yeasts and moulds in liquids or on surfaces.

Art. No.	Pack Qty.	Pack.
2956.1	20 unit(s)	box

## ROTI®DipSlide PCA/TSA-Letheen

For total germ count in liquids or on surfaces. Suitable for analysis of cosmetics and of preservative containing products or after disinfection. Contains two neutralizers.

Art. No.	Pack Qty.	Pack.
8715.1	20 unit(s)	box

## ROTI®DipSlide PCA/VRB

For total germ count and for detection of Coliforms in liquids or on surfaces.

Art. No.	Pack Qty.	Pack.
2925.1	20 unit(s)	box

## ROTI®DipSlide PCA

For total germ count in liquids or on surfaces.

Art. No.	Pack Qty.	Pack.
2927.1	20 unit(s)	box

## ROTI®DipSlide PCA/Legionella

For total germ count and for detection of legionells in solution or on surface.

-ROTI®DipSlide PCA/Legionella is suitable for the preliminary inspection of large-scale systems for heating drinking water, ventilation and air-conditioning systems, and water in evaporative cooling systems.

Art. No.	Pack Qty.	Pack.
2493.1	20 unit(s)	box

## ROTI®DipSlide PCA/VRBD

For total germ count and for detection of enterobacteria in liquids or on surfaces.

Art. No.	Pack Qty.	Pack.
2924.1	20 unit(s)	box

## ROTI®DipSlide CLED/MacCon

For detection of bacteria of the urinary tract and of enterobacteria/Coliforms in liquids or on surfaces.

Art. No.	Pack Qty.	Pack.
2957.1	20 unit(s)	box

## ROTI®DipSlide VRBD

For detection of enterobacteria in liquids or on surfaces.

Art. No.	Pack Qty.	Pack.
2954.1	20 unit(s)	box

Current prices at [www.carlroth.com](http://www.carlroth.com)

### Contact international:

Orders:  
Phone +49 721/5606 510 · Fax +49 721/5606 111  
[info@carlroth.com](mailto:info@carlroth.com) · [www.carlroth.com](http://www.carlroth.com)

Carl Roth GmbH + Co. KG  
Schoemperlenstr. 3-5 · D-76185 Karlsruhe

**Export prices might be higher.** All supplies and deliveries are subject to the terms and conditions of sale and delivery of Carl Roth GmbH + Co. KG, Karlsruhe

### Contact Schweiz:

Orders:  
+41 (0)61 / 712 11 60 · Fax orders: +41 (0)61 / 712 20 21  
[info@carlroth.ch](mailto:info@carlroth.ch) · [www.carlroth.ch](http://www.carlroth.ch)

ROTH AG  
Fabrikmattenweg 12 · 4144 Arlesheim

Net prices in CHF plus VAT. Orders over 200 CHF will not incur the flat shipping rate of 15 CHF. The General Terms and Conditions of Roth AG, Arlesheim, apply