



CLEAN ROOM
PRODUCT LINE

CLEAN ROOM PRODUCT LINE

The **CLEAN ROOM PRODUCT LINE** allows you to enter blender bags and sampling bags into your highly controlled environment easily and with complete confidence. Guaranteed sterile both inside and out thanks to a system of double pouch, it helps ensuring a clean environment and increase the efficiency of your operations.

The safe way to carry out tests and analysis in a clean room, without worrying about contamination.

Features and benefits

- Made of FDA compliant material
- Can accommodate solid, semi-solid and liquid samples
- Wide bag opening to facilitate insertion of sample
- Double sterility guaranteed
- Packaged with double pouch that allows to remove the outer bag before entering the clean room
- Suitable for use in the homogenizer
- RNase, DNase and pyrogen free




Inspired by your challenges

VISIT LABPLAS.COM FOR ADDITIONAL INFORMATION ABOUT OUR PRODUCTS | Tel: 1.450.649.7343

LABPLAS



CLEAN ROOM PRODUCT LINE BAG INSTRUCTION

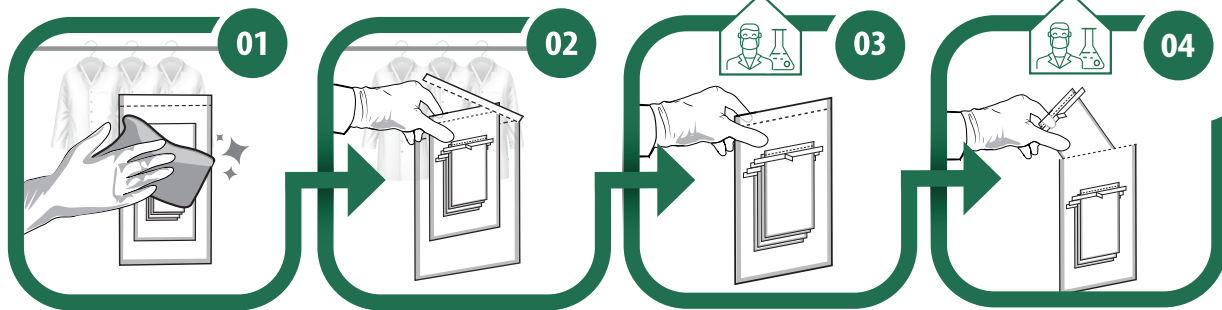


Once in the decontamination room, sanitize the first pouch

Remove the first pouch without contaminating the second pouch

Bring the sterile inner pouch into the clean room

Remove the inner pouch just before use

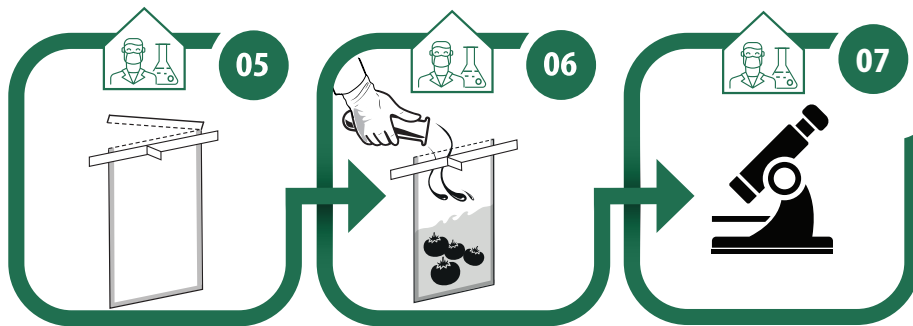

Tear off the top of the bag along the perforations

Aseptically insert the sample in the bag

Proceed according to the needs and specifications of the type of bag chosen for the analysis



Inspired by your challenges

LABPLAS

VISIT LABPLAS.COM FOR ADDITIONAL INFORMATION ABOUT OUR PRODUCTS | Tel: 1.450.649.7343

