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**Mercurisorb-ROTH®**

**Mercury binder**

Spilled mercury develops toxic, harmful vapors and should therefore be absorbed quickly and completely.

**Use:** For the quick absorption and easy absorption of spilled metallic mercury. Does not emit any more vapors after absorption.

**Available packs:**

**Emergency kits:**

Order no. : 9461.1 Emergency set in a plastic case 1 set  
Order no. : 9461.4 Emergency set in a box 1 set

An emergency set (in a plastic case or in a cardboard box) contains 100 g Mercurisorb-ROTH®, a small plastic scoop, card sheets, disposable Pasteur pipettes and an empty receptacle with a cap.

**Refill packs:**

Order no. : 9461.2 refill pack, small 250 g  
Order no. : 9461.3 refill pack, large 1 kg

**Contains: silver nitrate**

**Labelling:**

**Danger**

**Hazard statements**

H290 May be corrosive to metals  
H314 Causes severe skin burns and eye damage  
H410 Very toxic to aquatic life with long lasting effects

**Precautionary statements – prevention**

P273 Avoid release to the environment.  
P280 Wear protective gloves/protective clothing/eye protection/face protection.

**Precautionary statements - response**

P303 + P361 + P353 IF ON SKIN (or hair): Take off immediately all contaminated clothing. Rinse skin with water [or shower].  
P305 + P351 + P338 IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.

**Precautionary statements - disposal**

P501 Dispose of contents/container in accordance with local/regional/national/international regulations.

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EN

**Instructions for use:**

- Clear and cordon off the danger area, inform your superior.
- When spilled, fine mercury globules form, which can spread widely.
- Caution:** There is a high risk of spreading to other areas through shoes and objects.
- When cleaning up spilled mercury, wear protective goggles, protective gloves and, in the case of large quantities, respiratory protection.
- First push together larger mercury spheres using a suitable tool. Do not wipe or rub!
- Pick up as much as possible with the enclosed disposable pipette and fill it into the white container.
- Sprinkle Mercurisorb-ROTH® on very small or inaccessible mercury spheres in approximately the same weight.
- Mix Mercurisorb-ROTH® and mercury with the card sheets, avoid creating dust and keep moisture away (stains!).
- The absorption process can be clearly followed by the progressive gray coloration of the mixture. The process is over after a few minutes.
- Pick up with the enclosed scoop and pour into the plastic bottle.
- Label the receptacle and dispose of it for chemical disposal.
- Then ventilate the affected rooms generously.

**Annotation:**

The end product is a stable mercury compound that practically emits no more vapors. If stains have occurred due to the ingress of moisture, these can be removed with the special cleaner and stain remover Silvosol (order no. 0047). In the case of sensitive material, a resistance test should be carried out beforehand, if possible in an invisible area. It can be wiped or rinsed with water.

**Absorbance: 1:1,**

i.e. 100 mg Mercurisorb-ROTH® can bind 100 mg metallic mercury.

**Shelf life:** In the closed original packaging, at least 2 years from the date of production.

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